
CONGEN

SureTaq Hotstart Taq Polymerase

Art. No. F4005
100 rxn

User Manual

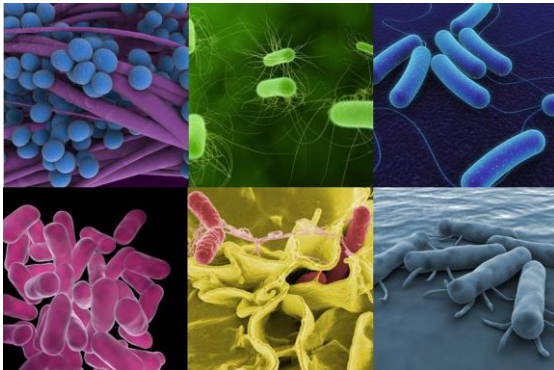January 2024

Inhalt

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1 General Information

1.1 Description

SureTaq Hotstart Taq Polymerase is an additional component for use with SureFast® / SureFood® real-time PCR Kits.

SureTaq Hotstart Taq Polymerase is not verified lot-specifically with SureFast® / SureFood® real-time PCR Kits.

1.2 Kit components and storage

| Kit Code | Reagent | Amount | Lid Color |
|----------|----------------|-----------|-----------|
| T | Taq Polymerase | 1 x 15 µl | Red |

Store the Taq Polymerase at **-20°C** and protected from light.

1.3 Additionally required equipment and materials

- DNA extraction kit (e.g. SureFast® PREP products / SureFood® PREP products)
- Real-time PCR methods (e.g. SureFast® products / SureFood® products)

2 Qualitative Analysis

2.1 Protocol

2.1.1 Preparation of the master-mix

Further information about the calculation of the total number of reactions needed (samples and control reactions) for the specific PCR assay is given in the manuals of the SureFast® / SureFood® real-time PCR products.

It is also recommended to prepare the master-mix with 10 % additional volume in order to compensate reagent loss. Allow the reagents to thaw, mix and centrifuge before opening and use.

Example for the calculation and preparation of 10 reactions:

| Components of the master-mix | Amount per reaction | 10 reactions (with 10% excess) |
|------------------------------|---------------------|--------------------------------|
| Reaction Mix | 19.9 µl | 218.9 µl |
| Taq Polymerase | 0.1 µl | 1.1 µl |
| Total volume | 20 µl | 220 µl |

Mix each master-mix well and centrifuge shortly before use.

3 Further Information

3.1 Technical Support

For further questions please send an e-mail to info@congen.de.