
CONGEN

SureFast® Escherichia coli Serotype I 4plex

Art. No. F5167
100 rxn

User Manual

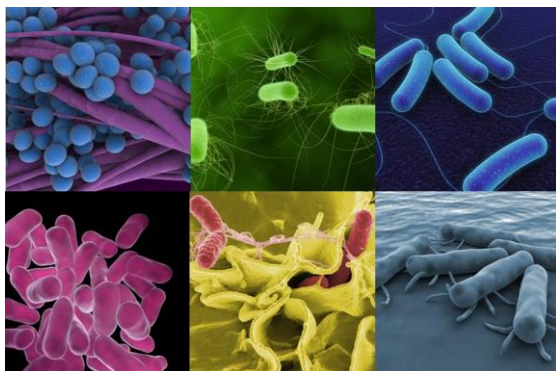June 2026

Inhalt

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Content

1	General Information	11
1.1	Description	11
1.2	Limit of Detection	11
1.3	DNA-preparation	12
1.4	Kit components and storage	12
1.5	Additionally required equipment and materials	12
1.6	Precautions for users	12
1.7	Setup.....	13
1.8	Detection channel Set-up	13
2	Qualitative Analysis	15
2.1	Protocol	15
2.1.1	Preparation of the master-mix	15
2.1.2	Preparation of the real-time PCR-mix	16
2.2	Interpretation of results	16
3	Validation	18
4	Limitations of the method	18
5	Further Information	18
5.1	Product Information	18
5.2	Technical Support	18

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1 General Information

1.1 Description

The SureFast® *Escherichia coli* Serotype I 4plex is a real-time PCR for the direct, qualitative detection and differentiation of specific *wzx* gene DNA sequences of *Escherichia coli* O121, O26 and O103 in enriched food and is intended to be used by all kinds of food testing laboratories.

SureFast® *Escherichia coli* Serotype I 4plex has been validated and certified in the context of an extraction using SureFast® STEC 4plex ONE (Art. No. F5265) or SureFast® PREP Bacteria (Art. No. F1021) by AOAC (PTM #052602).

Each reaction contains an internal amplification control (IAC). If the DNA contains PCR inhibiting substances, the signal of the amplification control will be affected or the amplification will be suppressed. Examples for PCR inhibiting substances are alcohols (e.g. ethanol, isopropanol), surfactants (e.g. CTAB, SDS, Triton X100) and salts (e.g. sodium chloride). In addition spices, herbs, algae, cocoa and further sample matrices might have PCR inhibiting effects.

The real-time PCR assay can be performed with commonly used real-time PCR instruments, equipped for detection of four fluorescence emissions at the channels FAM, VIC/HEX, ROX and Cy5 at the same time.

The external technical validation of instruments was performed on Bio-Rad CFX96 Deep Well DX, Bio-Rad CFX96 Opus Deepwell, and R-Biopharm RIDA®CYCLER.

The internal technical verification of instruments (not part of the PTM validation) was performed on Agilent Mx3005P, Agilent AriaDx, Applied Biosystems 7500, Roche cobas® z 480 Analyzer and Roche LightCycler® 480 Instrument II.

1.2 Limit of Detection

The SureFast® *Escherichia coli* Serotype I 4plex real-time PCR has a limit of detection of ≤ 5 DNA copies.

The assay limit of detection depends on sample matrix, processing grade, DNA preparation and DNA content.

The SureFast® PCR systems are very sensitive and therefore even a small amount of target DNA is sufficient for a successful analysis. The concentration of total DNA in the sample does not allow a conclusion on the quantity and quality of the target DNA.

1.3 DNA-preparation

For DNA-preparation the use of SureFast® PREP Bacteria (Art. No. F1021), Speed PREP (Art. No. F1054) or STEC 4plex ONE (Art. No. F5265) is recommended.

To assess the process of bacterial growth, it is suggested to compare the samples at the beginning and at the end of the culturing (bacterial growth at C_p difference > 3).

For cultivation the method of the DIN CEN ISO/TS 13136 is recommended.

1.4 Kit components and storage

Kit Code	Reagent	Amount	Lid Color
1	Reaction Mix	2 x 1050 µl	Yellow
2	Taq Polymerase	1 x 80 µl	Dark Red
3	Positive Control	1 x 200 µl	Light Blue

Store all reagents at -28 to -16°C and protected from light. The Taq Polymerase can be stored at +2 to +8°C for multiple uses on the same day.

Reagents should be used until the expiration date indicated on the label. Reagents can sustain up to 20 freeze/thaw cycles without influencing the assay performance.

Note: The Taq Polymerase may be in a frozen or unfrozen state. This does not affect the quality of the Taq Polymerase or the performance of the real-time PCR.

1.5 Additionally required equipment and materials

- DNA-Extraction kit
(e.g. SureFast® PREP Bacteria Art. No. F1021 / Speed PREP Art. No. F1054/ STEC 4plex ONE Art. No. F5265)
- real- time PCR instrument with four detection channels (510 nm, 580 nm, 610 nm and 660 nm)
- real-time PCR consumable (plates, tubes, foils, caps)
- pipettes with filter tips, for volumes from 0.7 µL – 1000 µL
- powder-free disposable gloves
- Vortex mixer
- Incubator
- Enrichment broth (i.e. modified Tryptone Soy Broth (mTSB) + Novobiocin)
- microcentrifuge with a rotor for the reaction tubes

1.6 Precautions for users

- Extraction, PCR preparation and the PCR run should be separated in different rooms to avoid cross-contaminations.
- This test must only be performed by laboratory personnel trained in molecular biology methods.
- Strictly follow the working instructions.
- When handling samples, wear disposable gloves. After finishing the test, wash your hands.
- Do not smoke, eat or drink in areas where samples or test reagents are being used.
- Food samples and enrichment cultures must be treated as potentially infectious as well as all reagents and materials being exposed to the samples and have to be handled and disposed according to the national safety regulations.
- Do not use the kit after the expiration date.
- All reagents and materials used have to be disposed properly after use. Please refer to the relevant national regulation for disposal.

SureFast® Escherichia coli Serotype I 4plex (100 rxn)

Art. No. F5167

June 2026

1.7 Setup

	Blockcycler ¹ & R-Biopharm RIDA®CYCLER ¹	Rotorcycler ² & LightCycler® 480 Instrument II ²
Initial Denaturation (HOLD)	1 min, 95°C	1 min, 95°C
Cycles	45	45
Denaturation	15 sec, 95°C	10 sec, 95°C
Annealing/Extension (CYCLE)	30 sec, 60°C	15 sec, 60°C
Temperature Transition Rate/ Ramp Rate	Maximum	Maximum

1.8 Detection channel Set-up

For additional details see cyclor operating instructions for real-time PCR device

Real-time PCR device	Detection	Detection channel	Quencher	Note
Agilent Mx3005P¹	<i>E. coli</i> O121	FAM	+	
	IAC	HEX	+	
	<i>E. coli</i> O26	ROX	+	
	<i>E. coli</i> O103	Cy5	+	
Agilent AriaDx / Mx¹	<i>E. coli</i> O121	FAM	+	
	IAC	HEX	+	
	<i>E. coli</i> O26	ROX	+	
	<i>E. coli</i> O103	Cy5	+	
Applied Biosystems 7500¹	<i>E. coli</i> O121	FAM	None	Check the passive reference option ROX is none.
	IAC	VIC	None	
	<i>E. coli</i> O26	ROX	None	
	<i>E. coli</i> O103	Cy5	None	
Bio-Rad CFX96/Dx/Opus (incl. Deep Well)¹	<i>E. coli</i> O121	FAM	+	Baseline Settings: - Baseline subtracted curve fit - Apply fluorescence drift correction
	IAC	VIC/HEX	+	
	<i>E. coli</i> O26	ROX	+	
	<i>E. coli</i> O103	Cy5	+	
R-Biopharm RIDA®CYCLER¹	<i>E. coli</i> O121	green	+	Ignore cycles before , if there is a significant deviation in the baseline at the start of the run. Please see page 45 of the cyclor operating instructions, section 12.1.2 Cycling analysis parameter.
	IAC	yellow	+	
	<i>E. coli</i> O26	orange	+	
	<i>E. coli</i> O103	red	+	

SureFast® Escherichia coli Serotype I 4plex
(100 rxn)

Art. No. F5167

June 2026

Real-time PCR device	Detection	Detection channel	Quencher	Note
Roche LightCycler® 480 II ²	<i>E. coli</i> O121	465-510	+	The SureCC Color Compensation Kit I (Art. No. F4009) is required.
	IAC	533-580	+	
	<i>E. coli</i> O26	533-610	+	
	<i>E. coli</i> O103	618-660	+	
Roche cobas® z 480 Analyzer ²	<i>E. coli</i> O121	465-510	+	The SureCC Color Compensation Kit I (Art. No. F4009) is required.
	IAC	540-580	+	
	<i>E. coli</i> O26	540-610	+	
	<i>E. coli</i> O103	610-670	+	

2 Qualitative Analysis

2.1 Protocol

2.1.1 Preparation of the master-mix

Calculate the total number of reactions needed (samples and control reactions) for the specific PCR assay.

Recommended control reactions for the specific PCR assay: negative control, extraction control and Positive Control.

For the analysis of enrichments additional controls are recommended: zero control and medium control.

The reaction mix contains an internal amplification control (Inhibition control) per reaction.

Reactions needed for the qualitative *E. coli* O121, O26 and O103 detection:

(e.g. when using samples without enrichment a.o. picked colonies or swabs - not part of the PTM validation* or after previous use of SureFast® STEC 4plex ONE, Art. No. F5265)

3 reactions for controls** (1x negative control, 1x extraction control, 1x Positive Control)

For each sample: at least 1 reaction for each sample DNA

Reactions needed for the qualitative *E. coli* O121, O26 and O103 detection in enrichments:

5 reactions for controls** (1x negative control, 1x extraction control, 1x Positive Control,
1x zero control, 1x medium control)

For each sample: at least 1 reaction for each sample DNA

It is also recommended to prepare the master-mix with 10% additional volume in order to compensate reagent loss. Allow the reagents to thaw, mix and centrifuge before opening and use.

Example for the calculation and preparation of 10 reactions:

Components of the master-mix	Amount per reaction	10 reactions (with 10% excess)
Reaction Mix	19.3 µl	212.3 µl
Taq Polymerase	0.7 µl	7.7 µl
Total volume	20 µl	220 µl

Mix each master-mix well and centrifuge shortly before use

* After DNA extraction with SureFast® PREP Bacteria (F1021), according to kit instructions.

**** Description of the controls**

- Negative control: only master-mix
- Extraction control: the extraction is performed without the sample – components from used Prep Kit
- Positive Control: master-mix and within the kit's provided Positive Control
- Zero control: master-mix and sample prior enrichment
- Medium control: only medium – no sample

2.1.2 Preparation of the real-time PCR-mix

- Pipette 20 µl of the master-mix into appropriate tubes/wells.
- Close the negative control.
- Pipette 5 µl of sample DNA into the designated tubes/wells and close them.
- Pipette 5 µl of Positive Control into the designated tubes/wells and close them.
- Centrifuge all tubes/plates shortly at low speed.
- Place tubes/plates into the real-time PCR instrument and start the run according to the setup.

2.2 Interpretation of results

The evaluation has to be made according to the usual analysis program recommended by the real-time PCR instrument manufacturer.

The control reactions have to show the correct results.

E. coli O121 DNA is detected in the FAM-channel, *E. coli* O26 DNA is detected in the ROX-channel and *E. coli* O103 DNA is detected in the Cy5-channel (see table). In the VIC/HEX-channel the internal amplification control (IAC) is detected.

A sample is stated **positive** for the respective parameter, if the sample DNA shows amplification in the respective channel. High amplicon concentrations can result in a weak or absent signal of the IAC.

A Cp value for the IAC is not needed to obtain a positive result of the Positive Control.

A sample is stated **negative** for the respective parameter, if the sample DNA shows no amplification in the respective channel and if the internal control (VIC/HEX-channel) of the sample is **positive**.

If the sample DNA in the VIC/HEX-Channel shows no amplification or sigmoidal amplification curve in the VIC/HEX channel, this indicates that inhibitors are present in the sample DNA that inhibit the PCR. Under these circumstances DNA isolation and purification of the sample need to be improved. When using a fast extraction it is recommended to centrifuge the lysate before the PCR or to repeat the extraction process. Alternatively the DNA can be diluted (recommendation 1:2 in PCR-water) and analysed again for inhibition. Please note that the dilution factor also affects the detection limit of the specific *E. coli* O121, O26 or O103 PCR assay.

Result in the respective channel				Interpretation
FAM channel <i>E. coli</i> O121	ROX channel <i>E. coli</i> O26	Cy5 channel <i>E. coli</i> O103	VIC/HEX channel IAC	
positive	negative	negative	positive/negative	<i>E. coli</i> O121 DNA detected
negative	positive	negative	positive/negative	<i>E. coli</i> O26 DNA detected
negative	negative	positive	positive/negative	<i>E. coli</i> O103 DNA detected
positive	negative	positive	positive/negative	<i>E. coli</i> O121 and <i>E. coli</i> O103 DNA detected
negative	negative	negative	positive	Negative, target DNA is not detected
negative	negative	negative	negative	invalid

Note: The results displayed in the table above represent merely an example. Additional combinations are also possible.

The following table shows the specification ranges of the kit controls

	Specification range	Internal amplification control (IAC)
negative control	negative	$28 \leq Cp \leq 36$
Positive Control (FAM – <i>Escherichia coli</i> O121)	$25 \leq Cp \leq 33$	not relevant
Positive Control (ROX – <i>Escherichia coli</i> O26)	$25 \leq Cp \leq 33$	not relevant
Positive Control (Cy5 – <i>Escherichia coli</i> O103)	$25 \leq Cp \leq 33$	not relevant

3 Validation

SureFast® Escherichia coli Serotype I 4plex was validated per the AOAC INTERNATIONAL Methods Committee Guidelines for Validation of Microbiological Methods for Food and Environmental Surfaces and the ISO 16140-2:2016 [Microbiology of the food chain -- Method validation -- Part 2: Protocol for the validation of alternative (proprietary) methods against a reference method].

The reference method for the matrix study was USDA Microbiological Laboratory Guidebook (MLG) 5C.04, "Detection, Isolation and Identification of Top Seven Shiga Toxin-Producing Escherichia coli (STEC) from Meat Products, Carcass, and Environmental Sponges" (USDA/MLG 5C.04) for raw meat matrices und US FDA Bacteriological Analytical Manual (BAM) "Chapter 4A: Diarrheagenic Escherichia coli" (FDA BAM Ch. 4a) for all other matrices.

The kit has been AOAC RI validated for the following matrices: raw beef trim, raw ground beef, baby spinach, sprouts, guacamole, wheat flower.

4 Limitations of the method

- The presence of PCR inhibitors may cause invalid results.
- Extremely low levels of target below the limit of detection (LoD) may be detected, but results may not be reproducible.
- In highly processed samples, the limit of detection may be shifted. Factors such as high pressures, mechanical stresses, chemical treatment, extreme temperatures and/or extreme pH values during manufacturing process – such as in canning production – can damage or degrade nucleic acids. This means that the sensitivity of the test kit may be reduced and not all original components may be detected.
- A positive test result does not necessary indicate the presence of viable organism. It is indicative for the presence of the target DNA (*Escherichia coli* O121, O26 and O103 DNA).

5 Further Information

5.1 Product Information

- Detailed information about setup of several real-time PCR devices (Download: www.congen.de/en/downloads)
- Product-related documents (Download: www.congen.de/en/eifu/)
- AOAC certificate (Download: www.congen.de/en/quality-standards/)
- Validation Report upon request

5.2 Technical Support

For further questions please send an e-mail to info@congen.de.